

AlteCap™ Septum capper

Quick start guide

Included in the box:

- AlteCap™ Septum capper
- Power supply
- Support adapters
- Quick start guide



The AlteCap Septum capper has been designed to work with most manufacturer's septum cap mats for sealing 2D barcoded with tubes in SBS racks, 24, 48, 96 well plates and deep well blocks.

Support adapters are included and are to be used if plates or racks require extra support during the capping cycle.

Unpacking instructions

1. Place the box on a flat surface and open the box
2. Remove the top foam support and the power supply
3. Remove the AlteCap™ Septum capper from the packing material and place the unit on to a level surface (e.g., lab bench).
4. Plug in the power supply and turn the AlteCap™ Septum capper on using the ON/OFF switch on the back of the unit

CAUTION: Please ensure that the unit is handled safely and that appropriate lifting methods are used. The unit weighs 12kg.

Using the AlteCap™ Septum capper with tubes in SBS rack

1. Place the septum cap mat on top of a rack of tubes
2. Ensure the septum caps align with the top of each tube
3. Place the rack of tubes and mat in the drawer



Septum cap mat

4. Press the **HIGH** button on the screen
5. Close the drawer. The drawer will lock once engaged
6. The drawer will automatically open once the capping cycle is completed
7. Remove the rack from the open drawer and peel back the mat. All tubes are now evenly capped



8. Follow the same instructions above when capping plates or deep well blocks

